

Software AUTOSIEVE



IDEAL FOR

- **AUTOMATIC SIEVE EVALUATION**
- **EASY MONITORING AND DOCUMENTATION OF YOUR SIEVING RESULTS**
- **EXTENSIVE PRESENTATION OF RESULTS**
- **AUTO-VALIDATION OF THE SIEVING PROCESS WITH THE VIBRATORY SIEVE SHAKER ANALYSETTE 3 PRO**



Software AUTOSIEVE

For automatic evaluation of sieving results

The extensive software AUTOSIEVE enables the automatic sieve evaluation with the Vibratory Sieve Shakers **ANALYSETTE 3 PRO** and **ANALYSETTE 3 SPARTAN** and the **Heavy Duty Analytical Sieve Shaker ANALYSETTE 18**. The software AUTOSIEVE is largely self-explanatory and enormously eases the weighing and the evaluation of your sieve analysis.

Simple, flexible, reliable.

AUTOSIEVE is easy to install and can be downloaded from www.fritsch-international.com/autosieve. All you have to do is to connect a laboratory analysis balance and then weigh the sieves before and after sieving using differential weighing. AUTOSIEVE will do the rest. The software enables the simple saving and retrieving of empty sieve weights, both of single sieves and complete sieve sets.

Especially clever by utilizing the Software AUTOSIEVE, the FRITSCH **Vibratory Sieve Shaker ANALYSETTE 3 PRO** can be controlled via a RS232-interface. All functions such as amplitude, sieving time, intermittent operation and type of sieving can be entered comfortably via the computer. The interface and AUTOSIEVE allow the monitoring of the sieving process via the online comparison of the set and actual amplitude and therefore an automatic validation of the sieving process.

FRITSCH Advantage: Test the Software AUTOSIEVE now, non-binding and free for charge for 90 days!



Download AUTOSIEVE free of charge now
www.fritsch-international.com/autosieve

Facts and Advantages:

- ⊗ Software AUTOSIEVE for all WINDOWS™
- ⊗ versions as of Windows 7 or higher
- ⊗ Auto-Validation of the sieving process with the FRITSCH Vibratory Sieve Shaker ANALYSETTE 3 PRO
- ⊗ Database for sieve sets and sieve tare weights
- ⊗ Control of various balances
- ⊗ Automatic weighing of individual sieves or complete sieve sets
- ⊗ Extensive presentation of the results: Presentation and printout as cumulative or frequency distribution and as graphs according to DIN ISO 9276-1
- ⊗ Pre-defined printout of results
- ⊗ Fixed value files for displaying quality relevant parameters
- ⊗ Display of reference curves
- ⊗ Subsequent input of already available sieving results
- ⊗ Different language modules
- ⊗ Recording and management of all parameters according to DIN 19683
- ⊗ Calculation of statistical values
- ⊗ Export of results to Excel™

Using Software AUTOSIEVE continuously

If you would like to continue to use the AUTOSIEVE programme after 90 days, just purchase an activation code: order no. 03.2902.00 from your merchant where you purchased your sieve shaker or contact us: www.fritsch-international.com/contact/.

After purchasing the software activation, you can install the AUTOSIEVE software on up to 3 Windows PCs.



ORDERING DATA

Order No. _____ Article _____

SOFTWARE AUTOSIEVE FOR WINDOWS

For controlling and automatic evaluation of the sieve analysis

- 03.2902.00 Software AUTOSIEVE-activation
for installation on up to 3 Windows PCs
(after expiration of the free of charge test phase)
incl. cable and device driver for controlling the
Vibratory Sieve Shaker ANALYSETTE 3 PRO
- 03.2600.00 Laboratory analysis balance up to 4.1 kg (± 0.01 g) with RS232
interface, incl. computer connection

Laboratory-analysis scale for the Heavy Duty Analytical Sieve
Shaker ANALYSETTE 18 on request

SYSTEM REQUIREMENTS – SOFTWARE AUTOSIEVE

AUTOSIEVE can be utilized with the Vibratory Sieve Shakers **ANALYSETTE 3 PRO** and **ANALYSETTE 3 SPARTAN** and with the **Heavy Duty Analytical Sieve Shaker ANALYSETTE 18**.

Software minimum requirements	Windows 7 or higher
Hardware minimum requirements	Standard WINDOWS™ PC with min. 2 GByte RAM